

# GXK-012


## 動物皿

### 取扱説明書

#### 適用機種

GX-Kシリーズ全機種

GF-Kシリーズ全機種

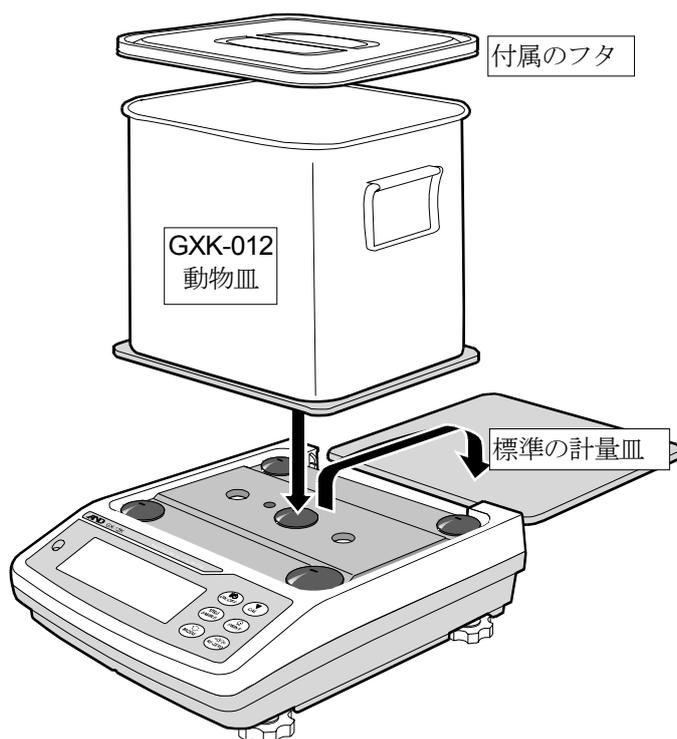

#### 使用方法

##### 準備

- 1 標準の計量皿を外し、GXK-012 動物皿に交換します。

##### 計量

- 2 **RE-ZERO** キーを押し、表示をゼロにします。
- 3 動物を入れ、計量します。

##### 注意


- 天びんのホールド機能（動物計量）を使用すると計量値の平均値を固定表示できます。
- 動物皿を使用した時はひょう量から約 1.5 kg 引いた値まで測定できます。
- 付属のフタは必要な場合に使用してください。
- キャリブレーションをする時は標準の計量皿を利用してください。
- 動物皿を載せた状態でキャリブレーションし、その後標準の計量皿を載せると **-E** 表示になります。標準の計量皿に交換する場合は標準の計量皿を載せた状態で再度キャリブレーションしてください。

# GXK-012

## Animal Weighing Bowl Instruction Manual

### Models

- GX-K series
- GF-K series

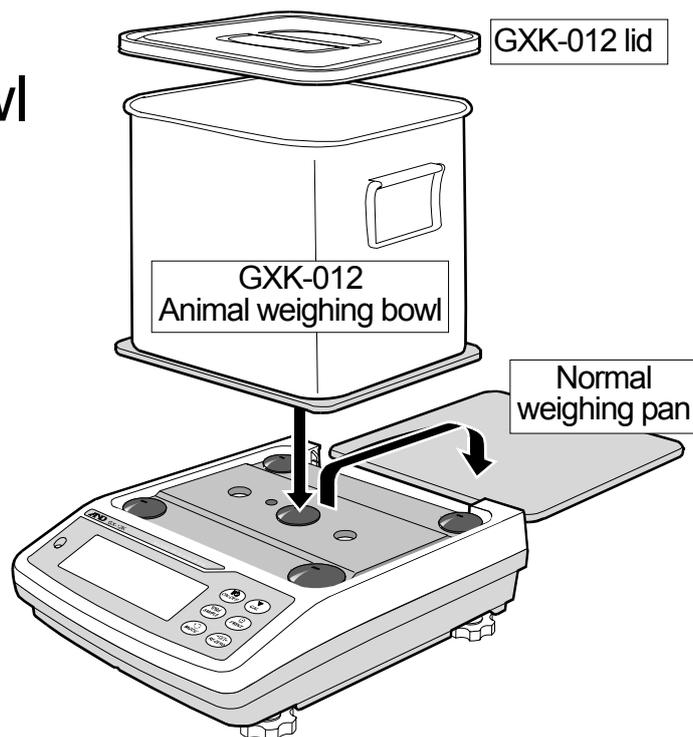

### Weighing

---

#### Setup

- 1 Replace the normal weighing pan with the animal weighing bowl (GXK-012).

#### Measurement

- 2 Press the **RE-ZERO** key to cancel the weight (tare).
- 3 Place a sample animal in the bowl to weigh it.

#### Caution

- When using the animal weighing mode (Hold function), the average weight of the animal can be displayed.
- When using this option, the weighing range that can be used is an approximately 1.5 kg less than the weighing capacity.
- If necessary, use the GXK-012 lid.
- When calibrating the balance, use the normal weighing pan.
- When calibrating the balance with the animal weighing bowl and replacing with the normal weighing pan, **-E** is displayed. If the normal weighing pan is used, re-calibrate the balance with it.